




Accurate, On-the-Spot Results

VICAM's AflaCheck® is a qualitative one-step test kit for the detection of aflatoxin. Aflatoxin is a Group 1 carcinogen proven to cause cancer in humans. In the United States and abroad, aflatoxin levels in food and grain are subject to strict regulations. The European Union has established regulations for aflatoxin in groundnuts; tree nuts, such as almonds and pistachios; dried fruits; spices; and cereals. The U.S. Food and Drug Administration has instituted maximum levels for aflatoxin in many commodities, including grains and groundnuts.

AflaCheck Test Strips provide fast, accurate, on-the-spot results for the presence of aflatoxin at detection levels of either 10 ppb or 20 ppb. AflaCheck offers the option of two testing procedures to meet your specific needs.

AflaCheck is easy to use and provides reliable results in as little as 3 minutes, works in the field or in the lab, and doesn't require the use of expensive equipment. Simply dip the test strip into the sample and read results. Used alone, or in conjunction with VICAM quantitative immunoaffinity columns, AflaCheck is an ideal first step for detecting aflatoxin contamination.

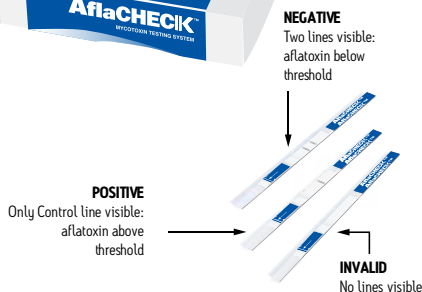

BENEFITS

Fast – Results in as little as 3 minutes*

Simple – No special training or equipment required

Accurate – Detects the presence of aflatoxin at levels of 10 ppb or 20 ppb

Convenient – Can be performed anywhere with a variety of samples

Durable – Long shelf life, no refrigeration required

Economical – Inexpensive first step in your testing protocol

Versatile – Choice of two cut off procedures; 10 ppb and 20 ppb

CULTIVATING SUCCESS THROUGH SCIENCE™:

1. **EXTRACT** sample with methanol and water.
2. **DILUTE** the extract with water.
3. **INSERT** test strip into diluted extract.
4. **INTERPRET** results visually.
 - Appearance of both test line and control line means aflatoxin level is below the test cutoff.
 - Appearance of the control line alone means aflatoxin level is above the test threshold.
 - No visible control line means test is invalid.

*excludes sample preparation and extraction



ORDERING INFORMATION

Cat. No.	Description
100000173	AflaCheck® Test Kit (25 Tests)**
100000175	AflaCheck Vial (contains 25 AflaCheck Test Strips)
600000812	Strip Test Pipettor (25 per bag)
600000813	Strip Test Dilution Tube (25 per bag)
600000827	40 mL Extraction Tube (25 per bag)
G4047	10 mL Graduated Cylinder
600000937	Graduated Cylinder, 25 mL
G4036	Digital Timer

Also Required; Must be Sourced Locally

ACS Reagent-Grade Methanol

Distilled Water

** Contains AflaCheck Test Strips, Strip Test Pipettors, Strip Test Dilution Tubes and Extraction Tubes for 25 Tests

Subject to change without notice.

©2013 Waters Corporation. Waters, The Science of What's Possible, VICAM, AflaCheck and Cultivating Success Through Science are trademarks of Waters Corporation.

The analytical methods presented in this data sheet have been researched and developed by VICAM to be used exclusively with AflaCheck products. These methods have been validated in the VICAM laboratories to perform to the specifications indicated in the AflaCheck procedures. The user assumes all risk in using AflaCheck procedures and products. VICAM makes no warranty of any kind, expressed or implied, other than that AflaCheck products conform to Waters' printed specification and quality control standards. Waters will, at its option, repair or replace any product, or part thereof, which proves to be defective in workmanship or material. VICAM's undertaking to repair or service such products is exclusive and is in lieu of all other warranties whether written, oral, expressed, or implied, including any implied warranty of merchantability or fitness for a particular purpose. VICAM shall have no liability for anticipated or lost profits or any loss, inconvenience or damage whether direct, indirect, incidental, consequential or otherwise, to person or property, or for strict liability or negligence arising from or in connection with the use of these assay procedures or AflaCheck products.

Orders:

1848 N Deffer Drive
Nixa, MO 65714 USA

Tel: +1.877.228.4244
+1.417.725.6588

Fax: +1.417.725.6102

orders@vicam.com

Headquarters:

34 Maple Street
Milford, MA 01757 USA

Tel: +1.800.338.4381
+1.508.482.4935

Fax: +1.508.482.4972

www.vicam.com